
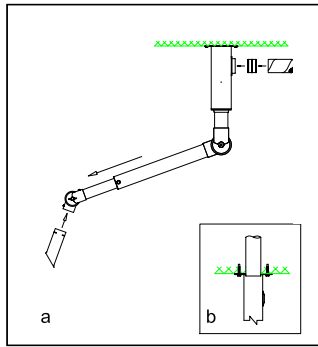
# CEILING MOUNTING – DECKENMONTAGE – LOFTMONTAGE

## UK – Teleskop – CM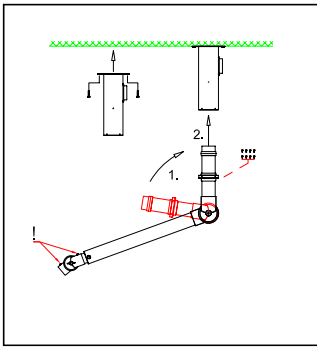



**Fig. 2**

- For side connection (fig. 2a) the main ventilation duct (125 mm dia.) can be connected to the outlet on the side of the column with a nipple – or directly to the outlet.  
For top connection the main ventilation duct is put through the ceiling and directly into the column. Make sure to seal the gap between the ceiling and duct with a filler.
- Loosen the fingerscrew located at the pipe reduction and pull the extraction arm to the length. Tighten the fingerscrew slightly again.
- Place the connection pipe of the accessory in the joint. Turn the slit on the connection pipe towards the small fingerscrew. When screwed in the fingerscrew will hold the accessory in place in the joint.
- Finally make sure the damper is open.
- The extraction arm is now ready for use.

- Considering the internal spring, it is important to turn the socket pipe in the direction of the arrow and that the fingerscrews (!) are faced forward.
- Place the socket pipe in the column. Turn the guide on the socket pipe until it falls into place in the column and fasten with the included screws.
- The extraction arm can now be turned 360° in the column.

Mount the column onto the deck or strut support frame. The arm operation will exert considerable strain and must be mounted on a solid surface.

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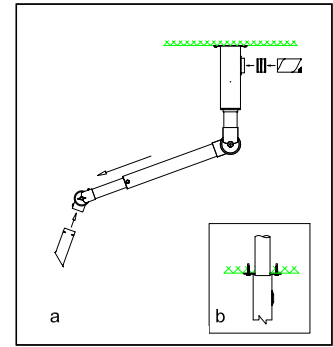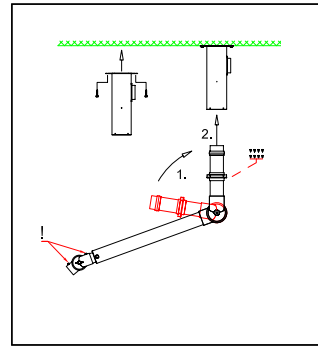
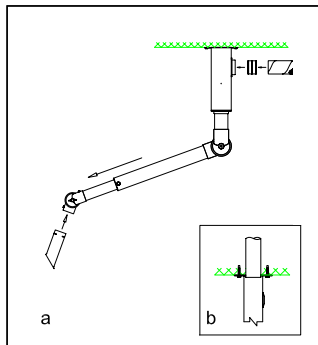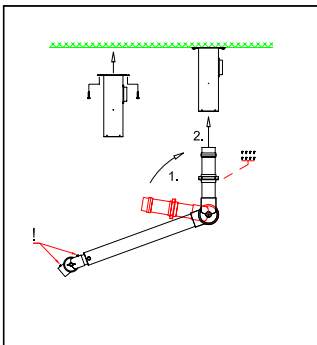
See MAINTENANCE AND CLEANING  
at [www.alsident.com](http://www.alsident.com) under "Technical Information"

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All rights reserved